



COSHH Assessment

Do you need to use a Fume Cupboard?

Are you using the correct type?

Training

Ensure you know how to use a Fume Cupboard safely and efficiently before starting your work.

Contact Area Safety Co-Ordinator



Check Air Flow

Panel on Fume Cupboard should read air flow $\geq 0.5\text{m/s}$

If Not Do Not Use

Sash Height

Sash should not be higher than 500mm when in use.

This improves Efficiency and reduces energy consumption.



Correct Set-Up

Minimise airflow disturbance

Minimise Equipment/Storage

Keep sill clear

Keep clean and tidy

Diversity Factor

Do you know what it means?

Do you know what to do if the alarm activates?



Emergency Response

Do you know what to do?

- Ventilation failure
- Chemical Spill
- Fire

Do you know who to report an incident to?

After Use

Dispose of waste correctly

Clean up any spillages

Remove unnecessary items

General Tidy

